



# Installation guide: FLUICS Print for eLabNext



**1** Unbox the printer, make sure the yellow box attached with the magnets to the top of the printer. Connect the cables in the following order:

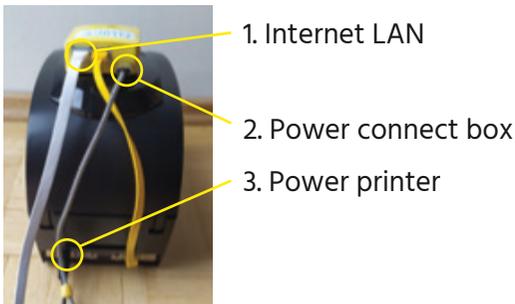


**2** Push the "Power" button to turn on the printer and wait a few minutes. **Check that all LEDs on the yellow box are on (except for the "status" LED that blinks every 3 seconds).**

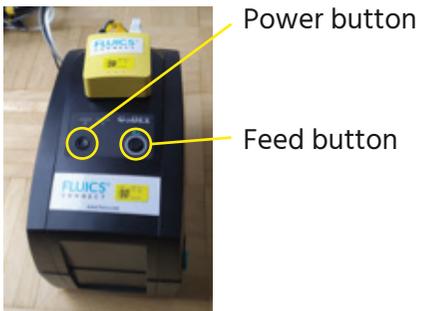


**3** Login to your eLabNext account install the FLUICS Print add-on from the Marketplace [elabjournal.com/members/marketplace](http://elabjournal.com/members/marketplace)



**4** Click "Configure" to enter the activation key which was sent to you by email. Didn't receive it? Write to [print@fluics.com](mailto:print@fluics.com)

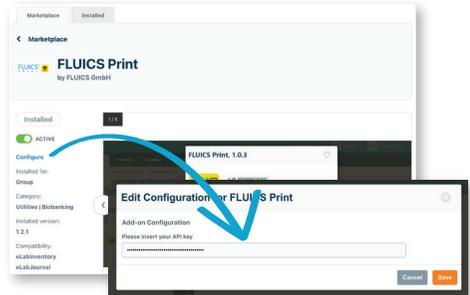


**5** Click on the "FLUICS" button to print labels for individual items

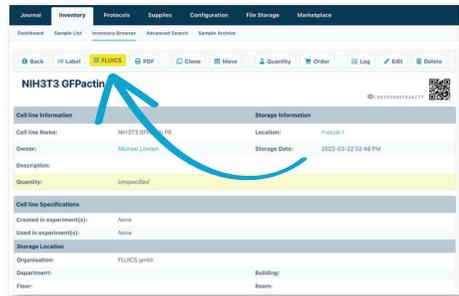

**6** Select label template and printer and click to print the label.



**7** Print labels for a selection of several items by clicking on the "Sample Actions" button.

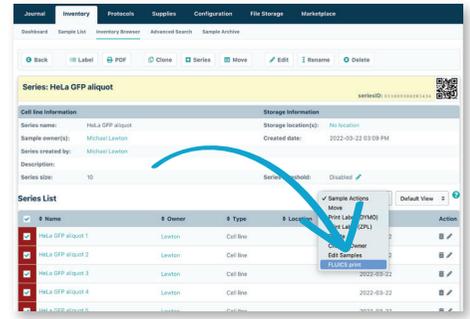

**Contact details:**

FLUICS GmbH  
 Fürstenstr. 15  
 80333 Munich  
 Germany

[connect@fluics.com](mailto:connect@fluics.com)  
 +49 89 71690877

Registered in  
 Munich HRB 246078